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PIPELINE  
TYPE

20

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SINGLE-USE LIQUID  
DISPENSING SYSTEM

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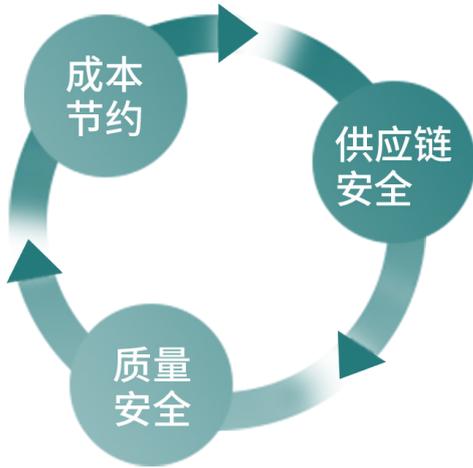
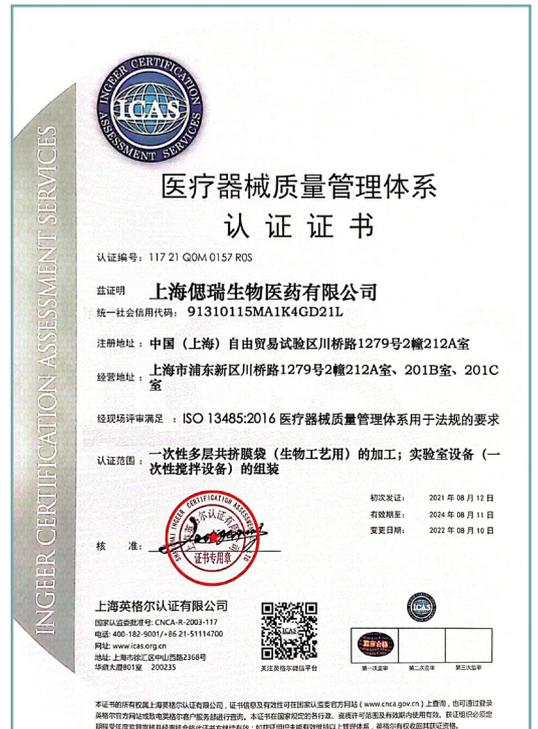
Shanghai Sirui Bio-Pharm Co., Ltd. is a high-tech enterprise engaged in the development, production and sales of single-use systems. Sirui is committed to building a high-end and high-quality domestic brand of single-use systems, and realizing the localization of single-use systems in various processes in the bio-pharmaceutical field.

Service philosophy: to provide customers with products and services of quality safety, supply chain security, and cost saving.

Quality and safety: Shanghai Sirui establishes and controls the system in accordance with the requirements of ISO9001 and ISO13485 standards. In order to meet the quality requirements of the client, our company manage the C-level production environment and the product realization process in accordance with GMP to ensure the stable and reliable quality of the products in each process and ensure the quality and safety of products.

Supply chain safety: Shanghai Sirui has effectively screened the suppliers according to the system requirements and the customized needs of the customers, and established a qualified supplier directory to ensure the supply of raw materials. Sirui conducts hierarchical audit management of suppliers and regular evaluation of the supply environment and quality control of raw materials to form a strategic partnership to ensure the safety of the product supply chain.

Cost saving: under the premise of ensuring quality, Shanghai Sirui gradually realizes the localization replacement of imported materials through development of raw materials. Sirui gradually reduces the cost of the industrial chain, and ultimately benefits the clinical benefit to the people."

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M

Operators perform and operate according to management documents.  
Work instruction for each department

M

Intact condition of equipment.  
Equipment management regulations  
Equipment validation

M

The quality of raw and auxiliary materials meets the  
production requirements.  
Procurement management rules  
Warehouse management regulations

M

Correct design, machining method and testing method.  
Product inspection procedure  
Production management procedure

E

Production environment in contact with the product, equipment hygiene,  
workshop tightness, workshop pressure, etc.  
Clean workshop environmental monitoring management regulations

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
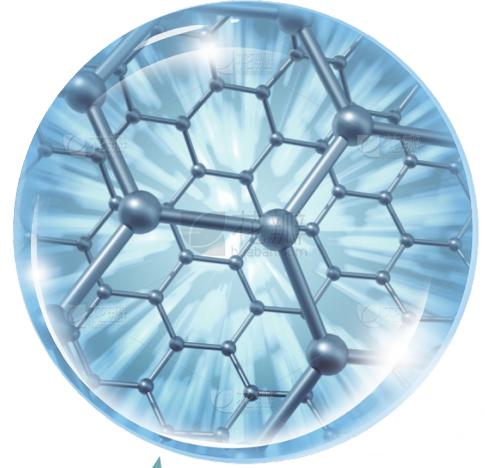
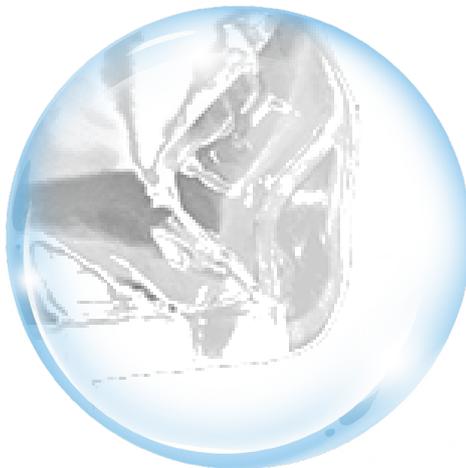
完整性测试  
Integrity Test

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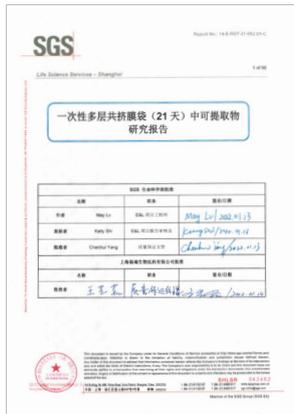Report of 21 days extractable study

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Report of the 70-day extractable study

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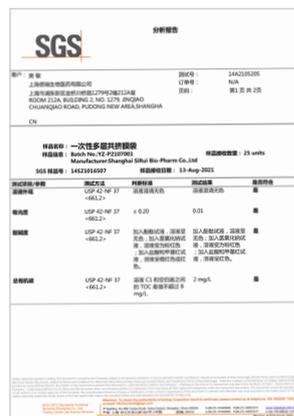
Chemical properties test (PH total organic carbon UV)

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
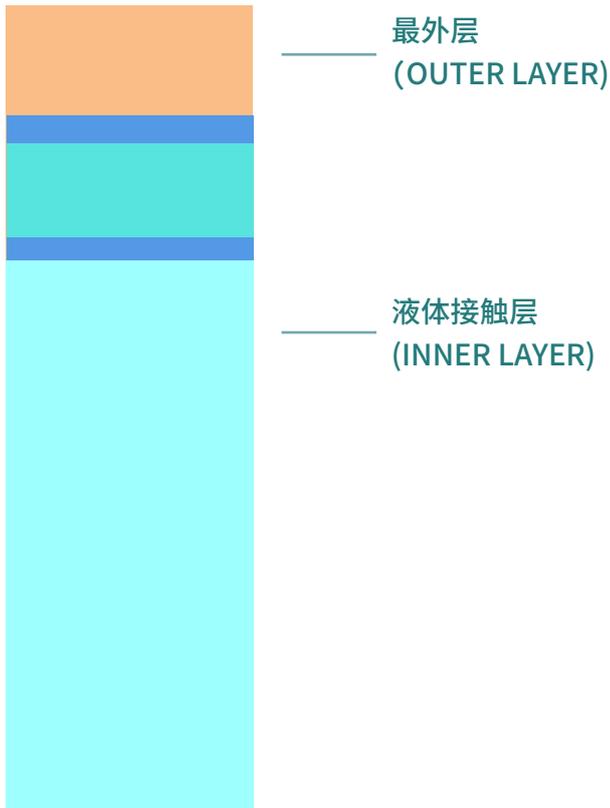
PE-tie-EVOH-tie-ULDPE（由外至内）厚度：0.325,mm

Disposable systems have extremely high requirements for materials, and the material of the bag body that directly contacts with the liquid on a large area is important. The 5-layer co-extruded film used in our bag making meets the relevant regulations.

The specific information is as follows:

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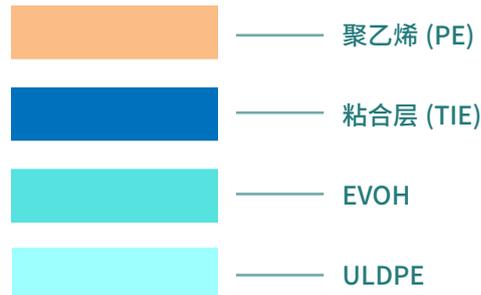**PE:** Excellent chemical compatibility and biocompatibility, good physical properties

**EVOH:** Super tear resistance, water and gas barrier layer, different permeability to various gases

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
## FILM REGULATORY

ISO10993-4:	Hemolysis
ISO10993-5:	Cytotoxicity
ISO10093-6:	Implantation test
ISO10993-10:	Irritation and Sensitization tests
ISO10993-11:	Acute Systemic Toxicity test
USP<85>:	Bacterial Endotoxins-LAL test
USP<88>:	Biological reactivity testing ,in vivo,class VI
USP<661>:	Plastic Containers European Pharmacopoeia tests, Ch. 3.1. 5
ADCF	

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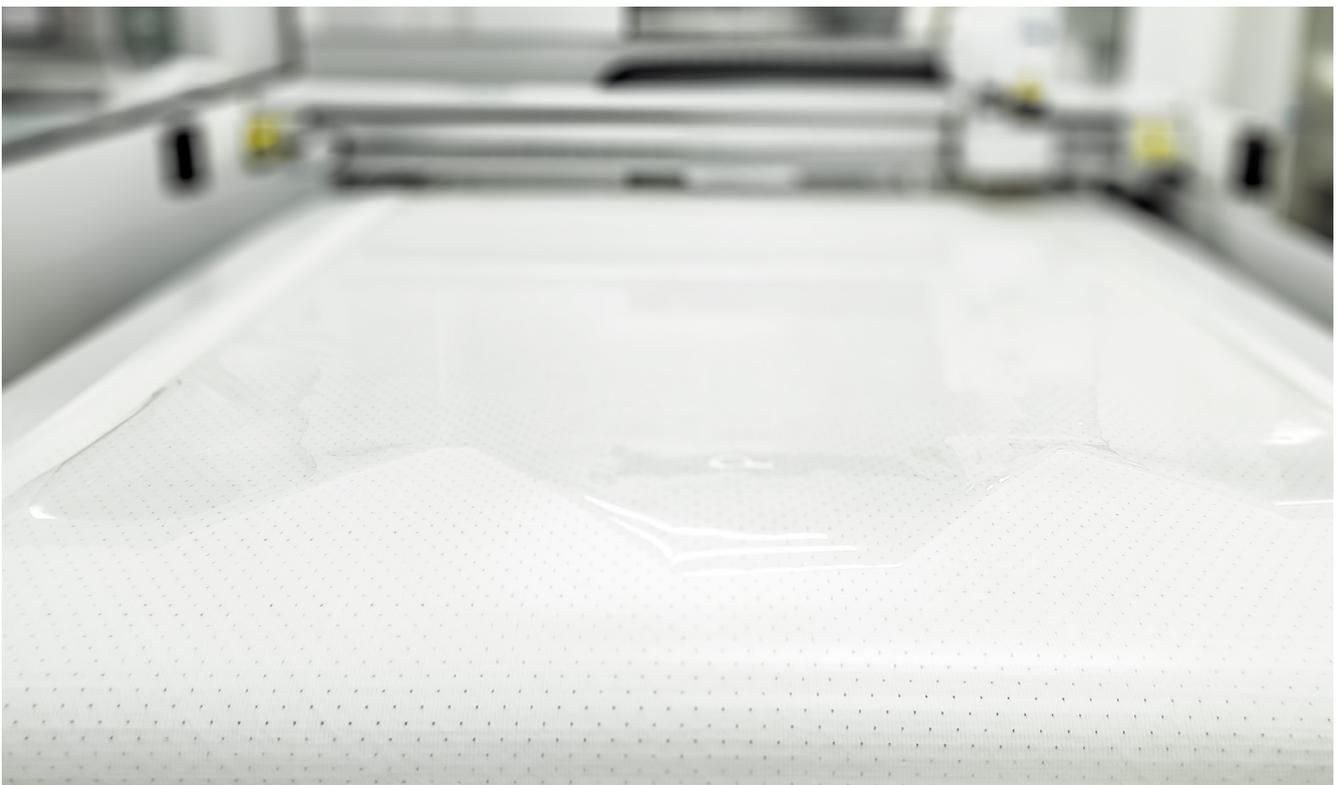The 2D Single-use Storage Bags produced by Shanghai SiRui Bio-pharm adopt multi-layer coextrusion film, which ensures the extremely low gas transmittance, good physical strength and extensive chemical compatibility and biocompatibility, so that it can be safely used for the storage and transfer of various biopharmaceutical liquids. Standard 2D Bags are available in sizes from 5mL to 50L, with special sizes to be customized. The shape design is divided into pipe type and boat type.

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## APPLICATION

- > Storage of media, feed, etc.
- > Purification buffer storage
- > Collection of components or samples
- > Multi-batch product or liquid-material combination
- > Storage during transport

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## FEATURES

- > Membrane Material with good physical strength, chemical compatibility and biocompatibility
- > High transparency membrane, easy to observe
- > Flexibly customized with various sizes and meet the needs of multiple process points

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## BAGS INFORMATION

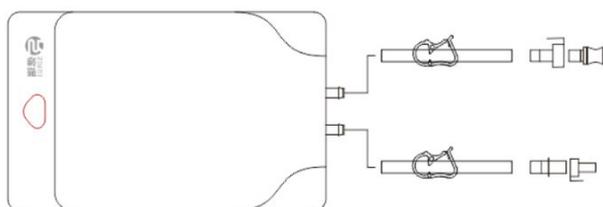
FILM	Multilayer film structure, liquid contact layer is ULDPE
TUBES	TPE, PVC, C-flex, etc., special specifications can be customized
CONNECTOR	Quick couplings, Luer sampling, filters, sterile connectors, Tube plugs, etc
TUBES NUMBER	2-3 Tubes for Standard products, and 1 Tube for 5mL
VOLUME	5mL-50L, with special specifications and sizes customized
STERILIZATION	Irradiation sterilization
PACKAGING	Two polyethylene outer layers
OPERATING TEMPERATURE	-80°C +45°C

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The pipeline type 2D liquid storage bag is connected by a number of 6.8\*8.8mm PE pipes to the bag body and the pipeline. The specification is 5ml-50L. The pipeline can be selected from various materials, apertures and lengths. The joints can be quick connector, Sterile connectors, Luer sampling, filters, pipe plugs, etc.

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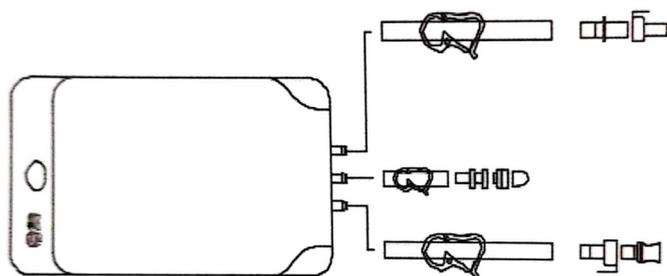
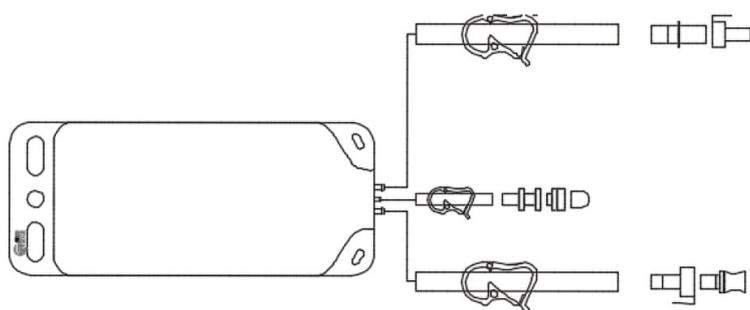
## ORDERING INFORMATION

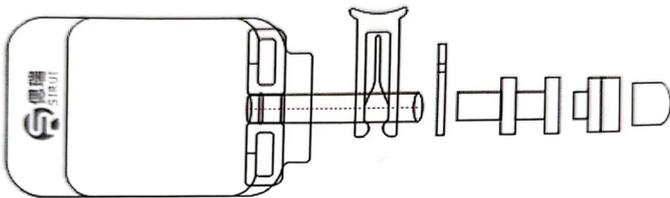
Article number	Volume	Pipeline	Packing specification
CX0005C	5L	3/8" X 5/8" TPE 30cm, MPC male+sealing cap, pinch clamp	30Pcs/Box
CX0010C	10L	3/8" X 5/8" TPE 30cm, MPC female+sealing cap, pinch clamp	15Pcs/Box
CX0020C	20L	1/4" X 3/8" TPE 10cm, Luer® Lock female + needle free	15Pcs/Box
CX0050C	50L	sampling port, dust cap, pinch clamp	15Pcs/Box

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The ship mouth 2D liquid storage bag is connected by the ship-shaped fittings to the bag body and the pipeline. The specification is 5ml-50L. The pipeline can be selected from various materials, apertures and lengths. The joints can be quick connector, sterile connectors, Luer sampling, filters, pipe plugs, etc.

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
### ORDERING INFORMATION

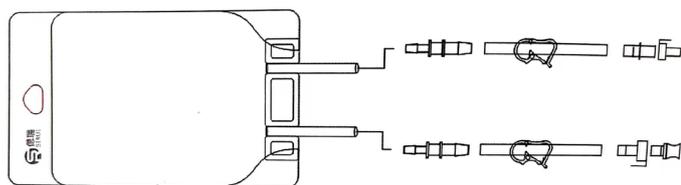Article number	Volume	Pipeline	Packing specification
2D005MA	5mL	6.8mmx8.8mm PE 5cm, Luer sampling, clip	200pcs/carton

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
## ORDERING INFORMATION

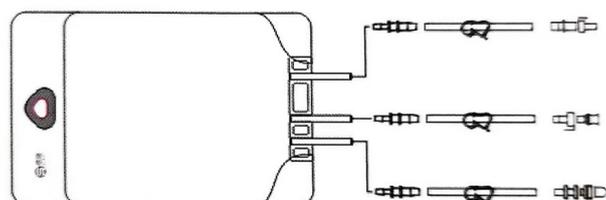Article number	Volume	Pipeline	Packing specification
2D050MB	50mL	1/8"x1/4" TPE 15cm, MPC male + sealing cap+pinch clamp	30pcs/carton
2D150MB	150mL	1/8"x1/4" TPE 15cm, MPC female + sealing cap+pinch clamp	30pcs/carton
2D250MB	250mL		30pcs/carton
2D500MB	500mL		30pcs/carton

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
## ORDERING INFORMATION

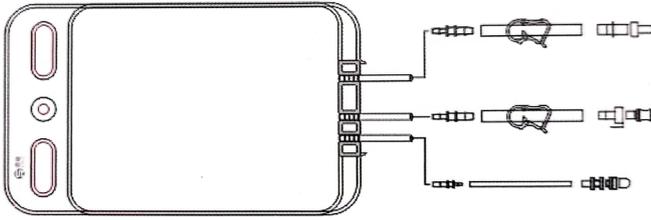Article number	Volume	Pipeline	Packing specification
2D0001C	1L	1/4" X 3/8" TPE 30cm, MPC male+sealing cap, pinch clamp	30pcs/carton
2D0002C	2L	1/4" X 3/8 " TPE 30cm, MPC female+sealing cap, pinch clamp	30pcs/carton
2D0003C	3L	1/8" X 1/4" TPE 10cm, Luer® Lock female + needle free sampling port, dust cap, pinch clamp	30pcs/carton

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
## ORDERING INFORMATION

Article number	Volume	Pipeline	Packing specification
2D0005C	5L	3/8" X 5/8" TPE 30cm, MPC male+sealing cap, pinch clamp	30pcs/carton
2D0010C	10L	3/8" X 5/8" TPE 30cm, MPC female+sealing cap, pinch clamp	15pcs/carton
2D0020C	20L	1/8" X 1/4" TPE 10cm, Luer® Lock female + needle free	15pcs/carton
2D0050C	50L	sampling port, dust cap, pinch clamp	15pcs/carton

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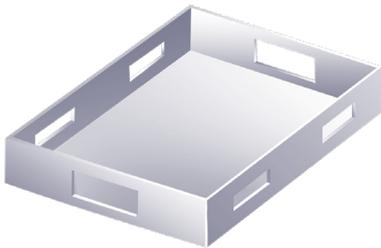### Plastic tray

Material	Polyethylene (PE)
Specifications	5L-50L (special specifications can be customized)
Features	<ul style="list-style-type: none"> <li>○ Ergonomic handle design</li> <li>○ Limit angle design, easy to maintain a stable stacking state</li> <li>○ The bottom plate is reserved for drainage pipe holes</li> </ul>
Application	2D storage bag assisted operation

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### Supporting stainless steel cart

Material	304 stainless steel (SS304)
Specification	Modular design, optional layer height/number of layers, easy for later expansion
Features	<ul style="list-style-type: none"> <li>○ Ergonomic handle design</li> <li>○ Side brackets can be used to hang liquid storage bags (≤ 10L)</li> <li>○ The baffle can be designed with a 3~5° inclination angle, which is conducive to liquid emptying</li> <li>○ Lightweight and strong, easy to transport</li> </ul>
Application	<ul style="list-style-type: none"> <li>○ 2D storage bag assisted operation</li> <li>○ 2D storage bag auxiliary operation</li> </ul>

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### Liquid storage bracket

Material	(SS304)
Specification	Modular design, optional layer height/number of layers, convenient Easy to expand later
Feature	<ul style="list-style-type: none"> <li>○ The baffle can be designed with a 3~5° inclination angle, which is conducive to liquid emptying</li> <li>○ Lightweight and strong, easy to transport</li> </ul>
Application	<ul style="list-style-type: none"> <li>○ 2D storage bag assisted operation</li> <li>○ Bulk intermediate product storage</li> </ul>

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Single-use liquid dispensing system

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The single-use liquid dispensing system is mainly composed of an upper mixing system, a disposable lining bag and a supporting container. It is suitable for uniform mixing and stirring of liquid-liquid and solid-liquid, which greatly improves the process efficiency of the process.

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### Upper mixing equipment

#### ● 304 stainless steel support:

to provide support for the entire platform during the mixing process

#### ● Imported motor and frequency converter:

accurate speed is convenient for verification of various mixing processes

#### ● Stainless steel mixing shaft and blades:

the installation and disassembly of the mixing blades is convenient and quick

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Matching container: optional plastic barrel or stainless steel barrel

Plastic bucket

- Linear low-density polyethylene (LOPE), resistant to common alcohol disinfectants
- The flared design is convenient for flanging and fixing of the lining bag
- The material is clean and uniform, the inner wall is smooth, and there is no damage to the soft bag body
- Maximum operating temperature is 60°C (no high temperature steam sterilization)
- Customize various specifications on demand
- Use with liner bag

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### Stainless steel barrel

- 304 stainless steel (SS304), resistant to wipe with common alcohol disinfectants
- The flared design is convenient for flanging and fixing of the lining bag
- The material is clean and uniform, the inner wall is smooth, and there is no damage to the soft bag body
- Customize various specifications on demand
- Use with liner bag

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### Supporting stainless steel transfer base

- Designed with matching container
- Main body: 304 stainless steel (SS304)
- Wheel surface: PU
- Pulleys: 4 silent universal wheels, all with locking function

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## PRODUCT INFORMATION

NUMBER OF INTERFACES	Customized
SPECIFICATION	19L-568L (special specifications can be customized)
OPERATING TEMPERATURE	-80°C ~+45°C
PACKING	Double independent packaging
STERILIZATION	γ-ray irradiation sterilization (25-40kGy)

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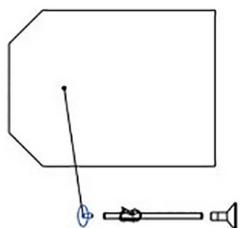
## FEATURES

- > Compared with closed bag type, it has higher economic efficiency
- > The open design is convenient for the rapid input of materials
- > The film material is highly transparent, which is helpful to assist the judgment of the process
- > Flexible customization of size, piping, joints
- > A variety of specifications, the volume can be customized 19L~568L

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## APPLICATIONS

- > Buffer preparation
- > Medium preparation
- > Intermediate product mixing

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Note: The connected pipes can be customized in various materials, diameters, and lengths, and the connectors can be quick connectors, bell mouths, etc.

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The 3D single-use storage bags produced by Shanghai SiRui Bio-pharm adopt multi-layer co-extruded film, which ensures the extremely low gas transmittance, good physical strength and extensive chemical compatibility and biocompatibility, so that it can be safely used for storage and transfer of various biopharmaceutical liquids. Standard 3D Bags are available in sizes from 50L to 3000L, with special sizes to be customized. SiRui 3D bioprocess bags can be used independently for the storage of buffers, media, intermediates or stock solutions; they can also be used in conjunction with a sterilization filter. Luer fittings, CPC quick connectors, tri-clamp sanitary fittings or sterile connectors are available for the inlet and outlet lines. We provide optional sampling methods, such as needleless samplers, Luer connectors, etc.

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## FEATURES

- > Membrane Material with good physical strength, chemical compatibility and biocompatibility
- > High transparency membrane, easy to observe
- > Flexibly customized with various sizes and meet the needs of multiple process points
- > Supporting a variety of tank or containers, compatible with multi-brand

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## BAGS INFORMATION

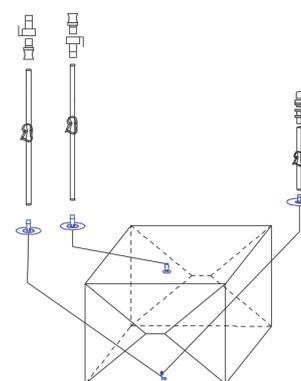
FILM	Multilayer film structure, liquid contact layer is ULDPE
TUBES	TPE, PVC, C-flex, etc., special specifications can be customized
CONNECTOR	Quick couplings, Luer sampling, filters, sterile connectors, Tube plugs, etc
TUBES NUMBER	1 Tube for inlet on the top, 2 Tubes for outlet and sample at the bottom
VOLUME	50L-3000L, with special specifications and sizes customized
STERILIZATION	Irradiation sterilization
PACKAGING	Two polyethylene outer layers
OPERATING TEMPERATURE	-80°C +45°C

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
## ORDERING INFORMATION

The 3D Single-use Storage Bags produced by Shanghai SiRui Bio-pharm Co., Ltd. are available in sizes from 50L to 3000L, with special sizes to be customized.

Part No.	3D0050C
Volume	50L
Top Port	3/8" X 5/8" Silicone 150cm, MPC female+sealing cap, pinch clamp
Bottom Port	3/8" X 5/8" Silicone 150cm, MPC male+sealing cap, pinch clamp
Sample Port	1/4" X 3/8" Silicone 20cm, Luer® Lock female + needle free sampling port, dust cap, pinch clamp

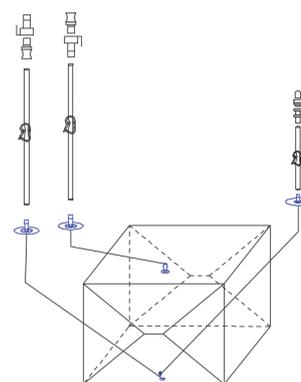
※Note: For 100 L, 200L volume 3D storage bags, Part No. are 3D0100C and 3D0200C.

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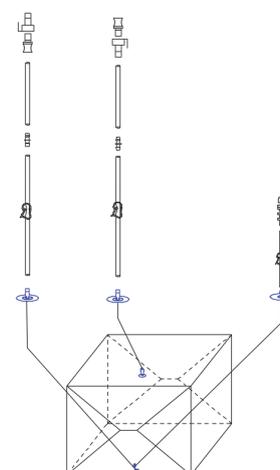
## ORDERING INFORMATION

<b>Part No.</b>	<b>3D0500C</b>
<b>Volume</b>	500L
<b>Top Port</b>	1/2"X3/4" Silicone 100cm+TPE 50cm+ MPX female + sealing cap, pinch clamp
<b>Bottom Port</b>	1/2"X3/4" Silicone 100cm+TPE 50cm+ MPX male + sealing cap, pinch clamp
<b>Sample Port</b>	1/4" X 3/8" Silicone 20cm, Luer® Lock female + needle free sampling port, dust cap, pinch clamp

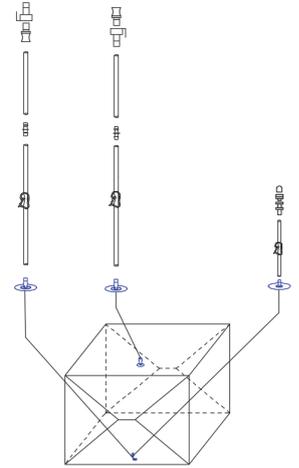※Note: For 1000 L, 2000L and 3000L volume 3D storage bags, Part No. are , 3D1000C, 3D1500C,3D2000C, 3D2500C and 3D2000C and 3D3000C.

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2D AUXILIARY PARTS FOR LIQUID STORAGE BAG

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### Foldable square storage box

<b>Material</b>	Polyethylene (PE)
<b>Specifications</b>	50L~1000L (special specifications can be customized)
<b>Feature</b>	<input type="checkbox"/> Strong and sturdy <input type="checkbox"/> Folding and stacking design, saving space and resources when not in use <input type="checkbox"/> Large size still has a comfortable height
<b>Application</b>	3D storage bag auxiliary operation and storage

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### Stainless steel square storage box

Material	304 stainless steel (SS304)
Specifications	50L~500L (Special specifications can be customized)
Feature	<input type="checkbox"/> Strong and sturdy <input type="checkbox"/> Detachable and stacked design, suitable for different solvents
Application	3D storage bag auxiliary operation and storage

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### Supporting stainless steel cart

Material	Main body: 304 stainless steel (SS304)
Specifications	Wheel surface: PU
Feature	to make according to the storage box size
Application	Used with storage box for 3D storage bag operation

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The same as 3D storage bags, 3D single-use mixing bags produced by Shanghai SiRui Bio-pharm adopt multi-layer co-extruded film, which ensures the extremely low gas transmittance, good physical strength and extensive chemical compatibility and biocompatibility, so that it can be safely used for storage and transfer of various biopharmaceutical liquids. Available in sizes from 50L to 3000L, mixing bags have a stirring paddle. They can be used in efficient mixing of buffers, media, intermediate or stock solutions; they can also be used in conjunction with a sterilization filter. Luer fittings, CPC quick connectors, tri-clamp sanitary fittings or sterile connectors are available for the inlet and outlet lines. We provide optional sampling methods, such as needleless samplers, Luer connectors, etc.

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## APPLICATION

- > Upstream culture medium, feed solid-liquid mixing
- > Buffer preparation and mixing for bioprocessing
- > Dissolution and mixing of hydrophilic/hydrophobic particles or powders
- > Suspension and mixing of bulk intermediates
- > Large volume terminal liquid dilution or concentration preparation

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## FEATURES

- > Membrane Material with good physical strength, chemical compatibility and biocompatibility
- > High transparency membrane, easy to observe
- > Suitable for bioprocess solid-liquid or liquid-liquid mixing
- > Supporting customization of special sizes or connectors, compatible with multiple brands

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## BAGS INFORMATION

FILM	Multilayer film structure, liquid contact layer is ULDPE
TUBES	TPE, PVC, C-flex, etc., special specifications can be customized
CONNECTOR	Quick couplings, Luer sampling, filters, sterile connectors, Tube plugs, etc
TUBES NUMBER	Regular includes a Feeding port, 1 Tube for inlet , 2 Tubes for outlet and sample
VOLUME	50L-3000L, with special specifications and sizes customized
STERILIZATION	Irradiation sterilization
PACKAGING	Two polyethylene outer layers
OPERATING TEMPERATURE	-80°C +45°C

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
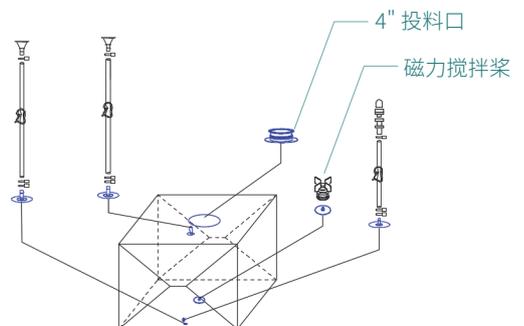
## ORDERING INFORMATION

The 3D Single-use Mixing Bags produced by Shanghai SiRui Bio-pharm Co., Ltd. are available in sizes from 50L to 3000L, with special sizes to be customized.

<b>Part No.</b>	<b>PD0050C</b>
<b>Volume</b>	50L
<b>Top Port</b>	3/8" X 5/8" Silicone 150cm, TC Sanitary fitting, pinch clamp
<b>Bottom Port</b>	3/8" X 5/8" Silicone 150cm, TC Sanitary fitting, pinch clamp
<b>Sample Port</b>	1/4" X 3/8" Silicone 20cm, Luer® Lock female + needle free sampling port, dust cap, pinch clamp

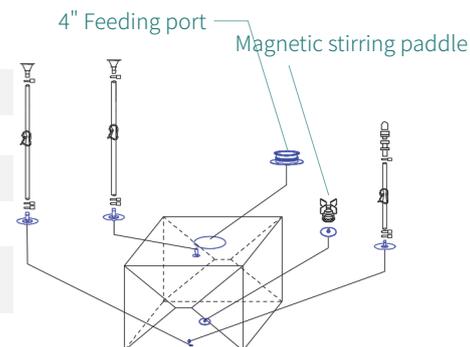※Note: For 100 L, 200L volume 3D Mixing bags, Part No. are PD0100C and PD0200C.


To order Mixing bags with temperature, pH or conductivity electrodes, please contact the technician or sales for customization.

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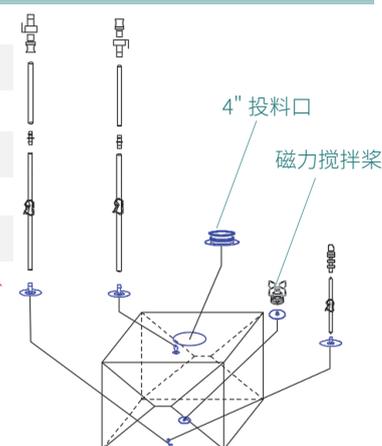
## ORDERING INFORMATION

<b>Part No.</b>	<b>PD0500C</b>
<b>Volume (L)</b>	500L
<b>Top Port</b>	1/2"X3/4" Silicone 100cm+TPE 50cm+ MPX female + sealing cap, pinch clamp
<b>Bottom Port</b>	1/2"X3/4" Silicone 100cm+TPE 50cm+ MPX male + sealing cap, pinch clamp
<b>Sample Port</b>	1/4" X 3/8" Silicone 20cm, Luer® Lock female + needle free sampling port, dust cap, pinch clamp

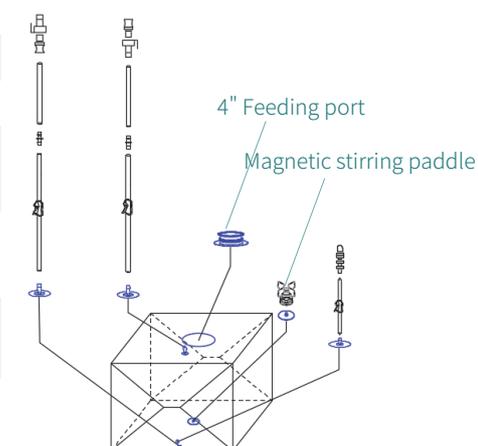※Note: For 1000 L, 2000L and 3000L volume 3D Mixing bags, Part No. are PD1000C, PD2000C and PD3000C.

To order Mixing bags with temperature, pH or conductivity electrodes, please contact the technician or sales for customization.

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The Single-use Magnetic Mixing System of Shanghai Sirui Bio-Pharm Co., Ltd. is driven by bottom magnetic force device and used to achieve powder-liquid or liquid-liquid preparation and mixing of various bio-pharmaceutical processes with matching 3D single-use mixing bags. Sirui Bio-Pharm's single-use mixing system can more efficiently and conveniently complete the preparation and mixing of media, feeding, additives, buffers, intermediates or concentrate. Compared with the traditional stainless steel mixing system, it can greatly save the manpower and material costs and time of SIP or CIP, save the cost of equipment cleaning and verification, and finally achieve efficient completion of the bio-process, and improve the turnover efficiency of the production line or equipment.

Sirui Bio-Pharm's single-use mixing system is mainly composed of a pre-assembled 3D single-use mixing bag with a built-in stirring paddle, an external support container, and mixing system controller with optional temperature control unit (TCU) module. Our mixing system can be configured with online weighing module and pH, conductivity or DO online monitoring modules, and data printing module. The standard mixing system volume is from 50L to 3000L, and offer special size and modules customization services, to cover preparation and mixing requirements of upstream and downstream bio-process, which can easily achieve stable amplified mixing effect.

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## APPLICATION

- > Buffer preparation
- > Media preparation
- > Intermediate products mixing
- > Preparation or concentrate of semi-products

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## FEATURES

- > Touch-screen operation interface, with perfect data integrity
- > Magnetically coupled stirring to achieve contactless stirring
- > Central control connection can be carried out to achieve central control integration
- > Pre-sterilized 3D single-use mixing bag, without SIP or SIP process
- > customizable, with selective modules to configure

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## ORDERING INFORMATION

The Single-use Magnetic Mixing System produced by Shanghai SiRui Bio-pharm Co., Ltd. ranges from 50L to 3000L, which are used in conjunction with the 3D Single-use Mixing Bags of SiRui. It is divided into basic and full models according to the different functional modules, and the actual functional modules can be specially customized according to the needs of users.

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## Basic Single-use Magnetic Mixing System

Volume(L)	Part No.	Length(mm)	Width(mm)	Height(mm)	Configuration Description
50	SR-CLJB-0050S	400	400	370	Magnetic stirring Module; Automatic control system (mixing software system, hardware electrical components); Operating system; Support tank (non-jacketed); Castors
100	SR-CLJB-0100S	500	500	450	
200	SR-CLJB-0200S	635	635	635	
500	SR-CLJB-0500S	835	835	835	
1000	SR-CLJB-1000S	1035	1035	1035	
1500	SR-CLJB-1500S	1275	1275	1005	
2000	SR-CLJB-2000S	2085	1120	1000	
2500	SR-CLJB-2500S	2160	1220	1020	
3000	SR-CLJB-3000S	2160	1220	1300	

For more configuration requirements, please contact the technical or sales.

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## Fully equipped Single-use Magnetic Mixing System

Volume(L)	Part No.	Length(mm)	Width(mm)	Height(mm)	Configuration Description
50	SR-CLJB-0050F	400	400	370	Magnetic stirring Module; Automatic control system (mixing software system, hardware electrical components); Operating system; Support tank (Jacketed); Castors; pH, DO and conductivity monitoring modules; Print module
100	SR-CLJB-0100F	500	500	450	
200	SR-CLJB-0200F	635	635	635	
500	SR-CLJB-0500F	835	835	835	
1000	SR-CLJB-1000F	1035	1035	1035	
1500	SR-CLJB-1500F	1275	1275	1005	
2000	SR-CLJB-2000F	2085	1120	1000	
2500	SR-CLJB-2500F	2160	1220	1020	
3000	SR-CLJB-3000F	2160	1220	1300	

For more configuration requirements, please contact the technical or sales.

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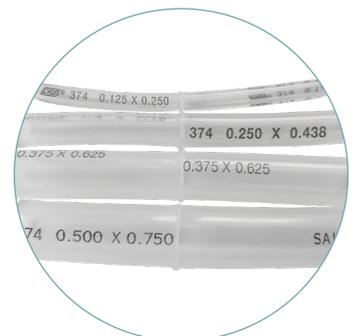
Tube-Welder I 是一种全自动装置，用于无菌焊接热塑管。  
The Tube-Welder I is a fully automated device for connecting thermoplastic tubing in a sterile welding operation.

该全自动装置可无菌连接外径 1/4 至 3/4 英寸的热塑管。  
The fully automated device allows sterile connection of tubing with outer diameter from 1/4" to 3/4".

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## FEATURES OF TUBE-WELDER

Easy to use. Reset operation can be executed even the upper cover is open. With one key reset and one key welding. Preinstalled settings for C-Flex®, AdvantaPure AdvantaFlex®, Pharmed® and Sanipure® tubing. True thermal weld provides exceptional strength. Aseptic connection without a laminar flow hood. Compact design, portable. Easy to validate, with validation equipment. One machine can handle tubes with different diameters.

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## Technical Specification

Tubing size	(ID X OD) 1/8" x 1/4"; 1/4" x 7/16"; 3/8" x 5/8"; 1/2" x 3/4" ect.
Tubing types	C-FLEX、PharMed BPT、Advantaflex ect. TPE Tubing
Welding cycle times	1-2 min (Depending on the tubing size and brand)
Dimensions (L*W*H)	260mm*248mm*248mm
Weight	9 kg
Material	Anodized aluminum
Positioning accuracy of three-axis motion	5 μm
Dimensions of blades	83mm*25mm*0.4mm
Use environment	Room temperature, 35-65% RH
Cooling mode	Forced air cooling
Power requirements	100-240V, 47-63 Hz
Rated power	288 W

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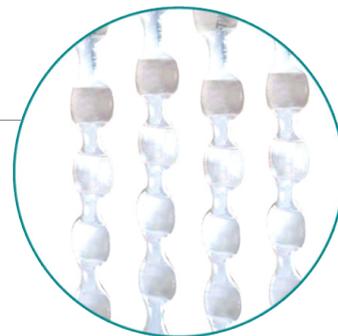
Tube-Sealer 是一种全自动装置，用于永久性、无泄漏密封热塑管。  
Tube-Sealer is a fully automated device, designed to produce permanent and consistent leak free seals on thermoplastic tubing.

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## TECHNICAL SPECIFICATION

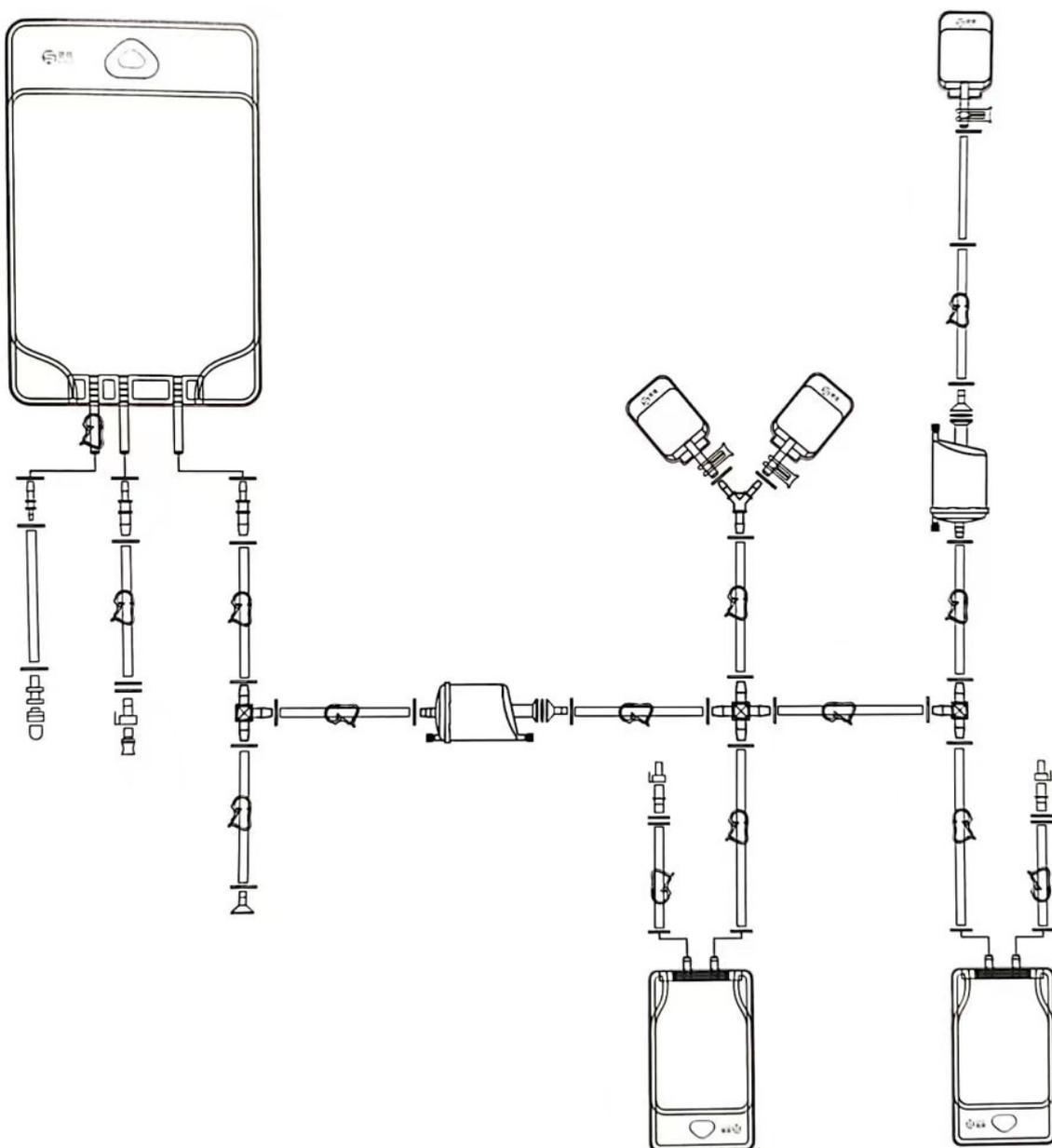
Tube size	Type I can handle tubing with max OD of 7/8", type II can handle tubing with max OD of 1.0") Wall thickness 1/16"-1/8" (1.6-3.2 mm)
Tube types	C-FLEX、PharMed BPT、Advantaflex TPE
Sealing cycle time	Depending on the tubing size and brand, sealing cycle of Tube sealer I is less than 4 min and sealing cycle of Tube sealer II is less than 5 min)
Dimensions (L*W*H)	210mm*165mm*135mm
Weight	About 3kg
Material and Surface Treatment	Aviation aluminum, anodized, it can be sterilized with 75% alcohol.
Use environment	Environment temperature: 15~35°C, Relative humidity 35-65%
Cooling mode	Forced air cooling
Temperature sensor	Pt temperature sensor
Program select mode	Select the sealing program by press the push button
Permissible pressure range	
liquid in tube	During sealing process: -1~+0.5 bar., After sealing process: -1~+1.0 bar, but not higher than specification of tube manufacturer. Tube has to be cooled off to room temperature.
Power requirements	100-240V AC, 47-63 Hz
Rated power	75W

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In practice, standard products may not be able to meet different usage requirements, so we have launched a convenient customized assembly and connection service. Commonly used accessories include various hoses, joints, sterile connectors, filters, liquid storage bags, etc.

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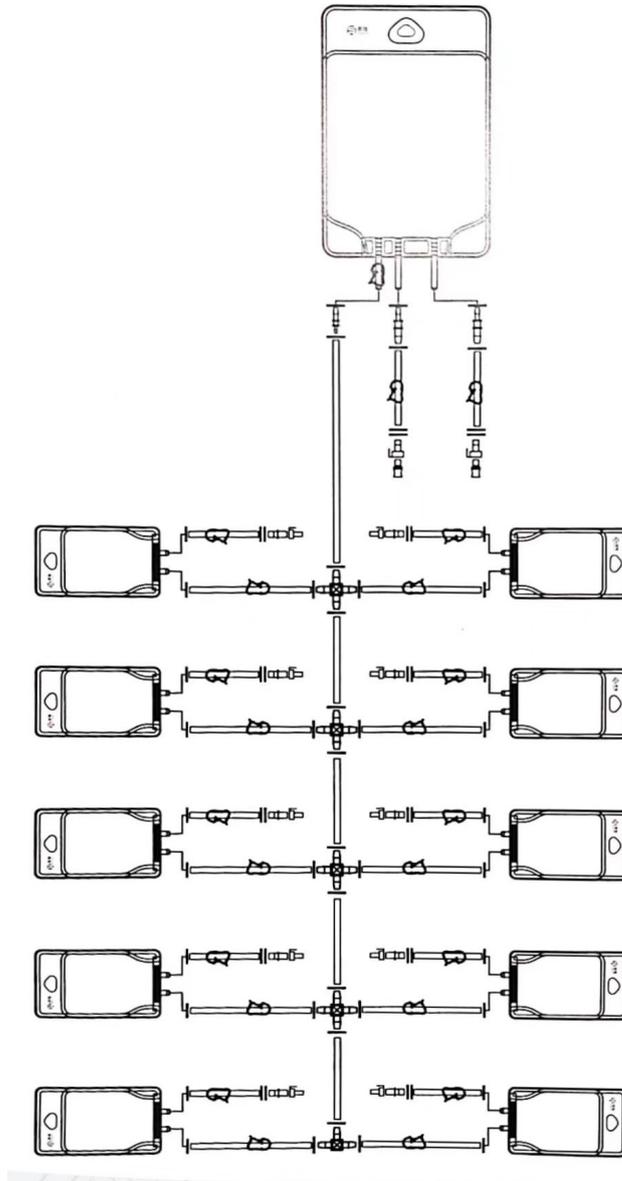
Each individual pipeline can be equipped with a sterile connector, and a sterile dis-connector; the pipeline can be quickly put into use after pre-irradiation and sterilization.

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